



TOC – Preparing your instrument for shutdown

1. Run a series of 3 DI water blanks through the instrument to flush it out.
2. Shutdown instrument through the software.
 - PC Controlled instruments:

Click  Shutdown

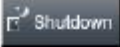

The Shutdown dialog box is displayed.

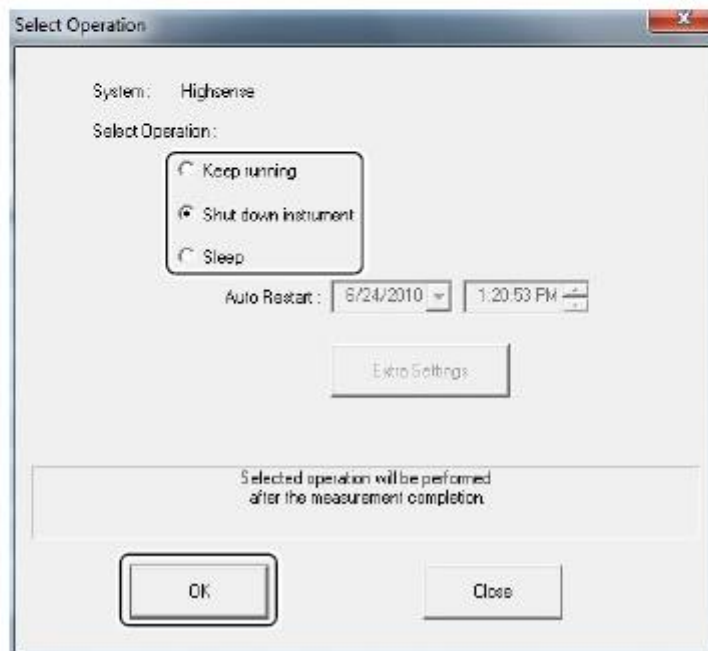

Push 'OK'. The instrument will start the shutdown procedure and after 30 minutes will completely shut itself off.

Use the main power button on the right hand side of the instrument to turn off power to the instrument.


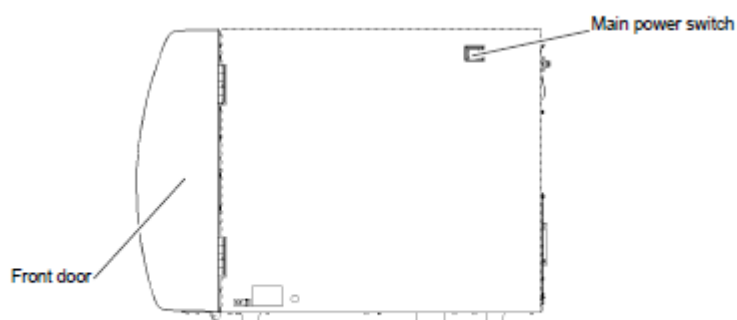


Figure 2.2 Right Side View

- Standalone Instruments:


Press the F1 [Standby Option] key in the "Initial Display".
The "Standby Option - Sleep" screen is displayed.

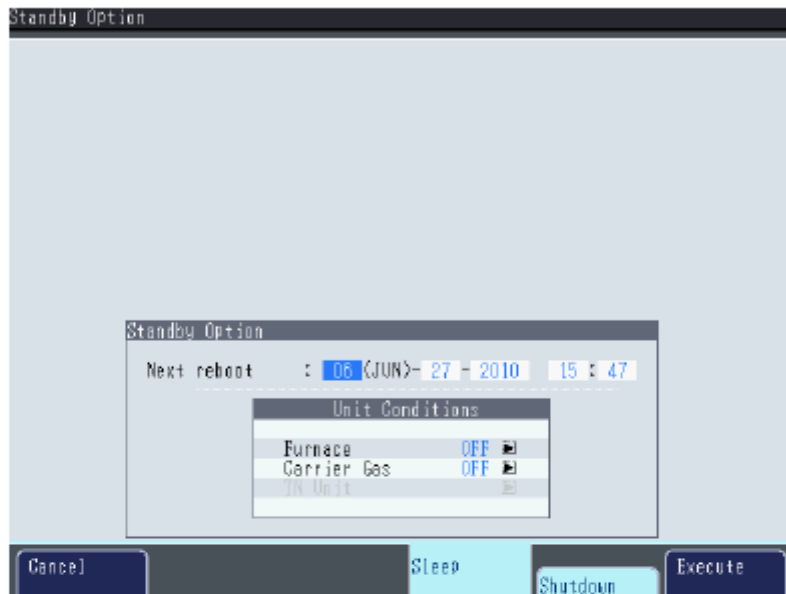

Press the F5 [Power OFF] key.
The message "Automatic Shutdown" is displayed.



Press the F6 [Execute] key.
"Do you want to stop this process?" is displayed.

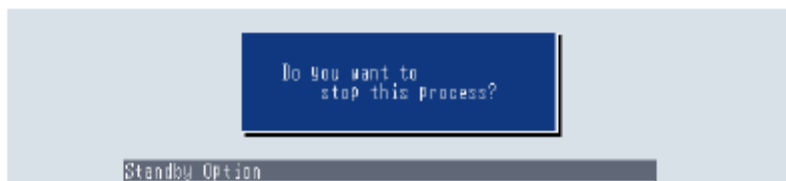

Press the F6 (Yes) key to confirm the shutdown. After 30 minutes the instrument will completely shut itself off. Use the main power button on the right hand side of the instrument to turn off power to the instrument.

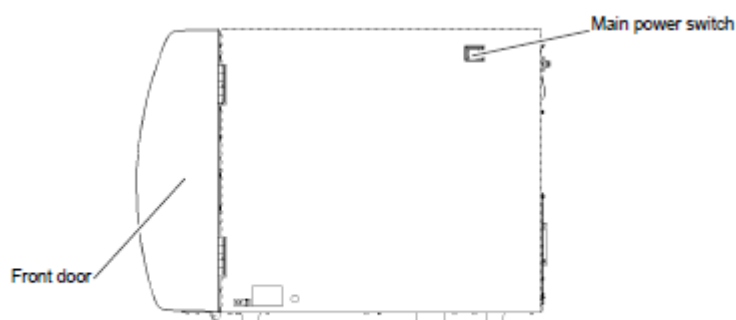


Figure 2.2 Right Side View

3. Close the Main carrier gas valve at the source.